

CALIBRATION INSTRUCTIONS
FOR NO. 2159 THERMOSTAT
MODEL 25X ALL-AMERICAN
PRESSURE STERILIZER

START STERILIZER AS IF YOU WERE GOING TO RUN A CYCLE.

IF UNIT SHUTS OFF AT TOO LOW OF TEMPERATURE OR PRESSURE, FOLLOW THESE INSTRUCTIONS:

TURN THERMOSTAT SHAFT COUNTER-CLOCKWISE. WHEN THE UNIT IS AT DESIRED TEMPERATURE OR PRESSURE, TURN SHAFT CLOCKWISE UNTIL LIGHT GOES OUT (DO NOT TURN ANY FURTHER). TIGHTEN THE KNOB ONTO THE SHAFT WITH POINTER IN THE UPRIGHT POSITION.

IF UNIT RUNS TOO LONG ON HIGH, FOLLOW THESE INSTRUCTIONS:

WHEN UNIT IS AT DESIRED TEMPERATURE OR PRESSURE, TURN SHAFT CLOCKWISE UNTIL LIGHT GOES OUT AND TIGHTEN KNOB IN UPRIGHT POSITION.