

# Macroscope Portable Microscope



*"a world of learning"*

This hand-held microscope is ideal for obtaining a close-up view of all kinds of subjects including biological specimens, textiles, coins, stamps and minerals. The clear acrylic base holds the microscope above the subject to allow fine focus and measurement using the scale built into the eyepiece lens. Being light-weight, portable and without the need for batteries, the Macroscope can be used almost anywhere to reveal the fine details of subjects that would otherwise be very difficult to see. A padded carry-case is included.

The top part of the Macroscope can be removed as used as a telescope, leaving the lower part as a magnifier.



## Instructions:

Position the subject below the acrylic base of the Macroscope and bring the measuring scale into focus by rotating the Upper Focus Ring. Bring the subject into focus by rotating the Lower Focus Ring.

Separate the components of the Macroscope by holding the Detachment Ring and turning the upper body anti-clockwise. The lower part can be used as a small magnifier and the upper part can be used as telescope.

## Specifications:

Product Code:	M3.20
Magnification:	45X
Working Distance:	42mm
Field of View:	4.8mm
Weight:	290g



Macroscope after detachment showing magnifier on left and telescope on right.

**Southern Biological - A Division of Cogitamus Pty Ltd**

ACN 097 468 954 ABN 25 838 926 933  
19-21 Worrell Street or PO Box 57 Nunawading Victoria 3131 Australia  
Phone: +61 3 9877-4597 Fax: +61 3 9894-2309  
Email: [sales@southernbiological.com](mailto:sales@southernbiological.com) Web: [www.southernbiological.com](http://www.southernbiological.com)